


Installing the Cobolt USB Laser Device Driver on a Windows 7 or Windows XP Machine


The USB configured controllers requires a driver for communication with a computer. The USB driver must be installed and can be downloaded from the Cobolt website (www.cobolt.se). When installed, a virtual COM port will be created to communicate with the Controller.

To install the USB driver in Windows 7 follow these instructions:


1. Go to the **Control Panel** and choose **Hardware and Sound**.

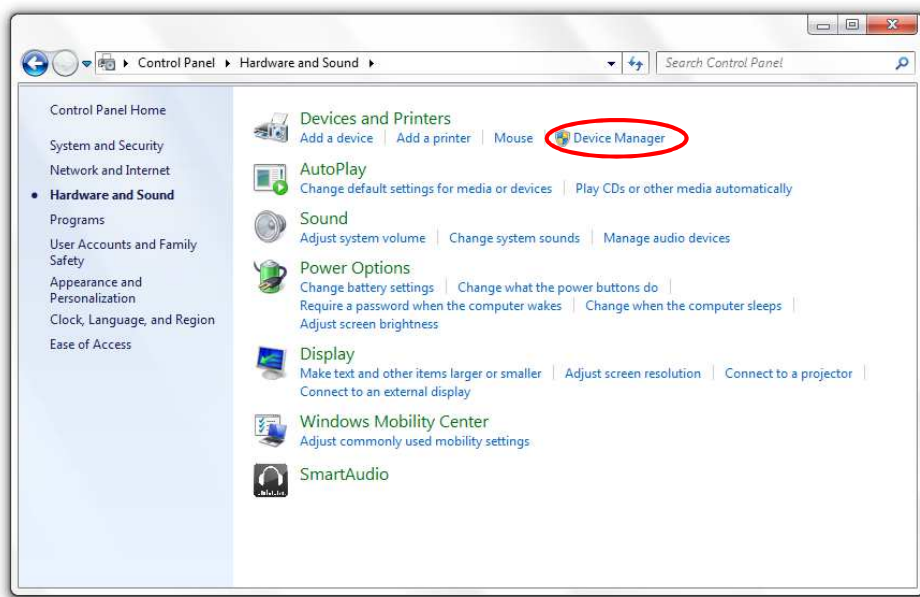


2. Under the **Devices and Printers** section, choose **Device Manager**.
3. Under **Other devices**, find the device called "**Cobolt Laser ...**". Right-click it and chose **Update Driver Software**.

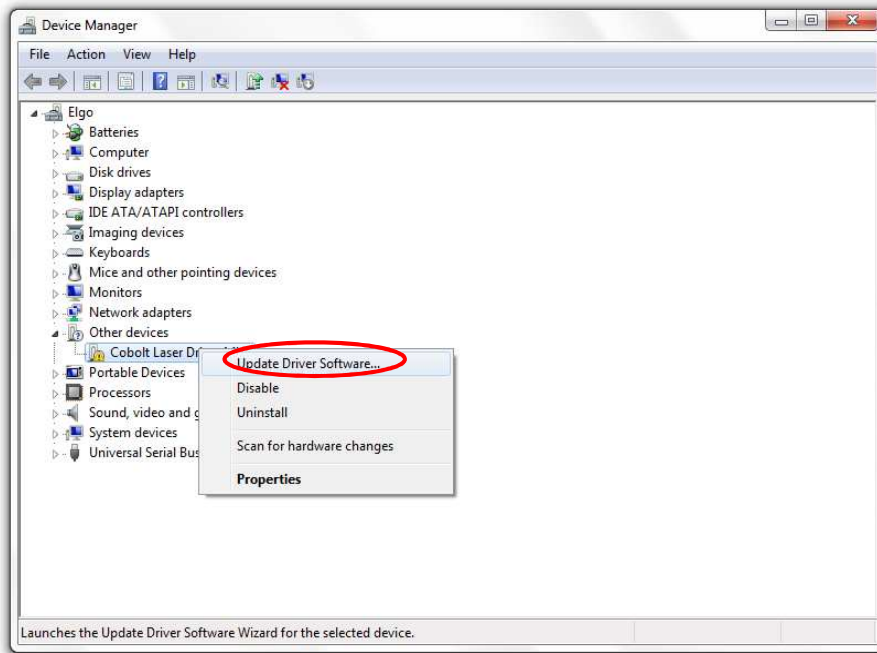

4. On the next screen chose the **Browse my computer for driver software** option.

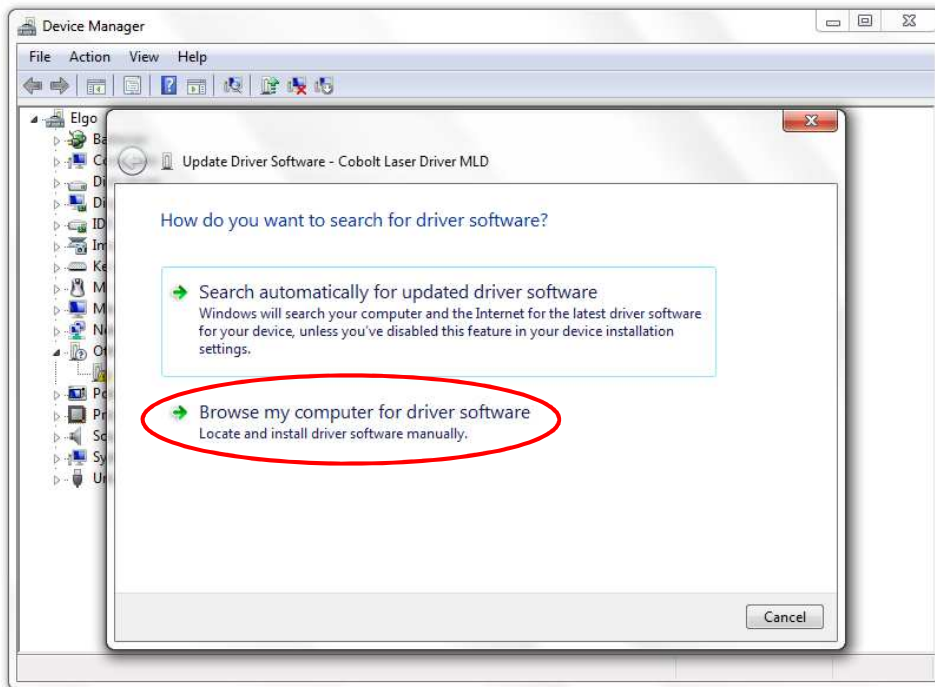

5. Click **browse**, and find folder on your computer where the USB driver is stored.

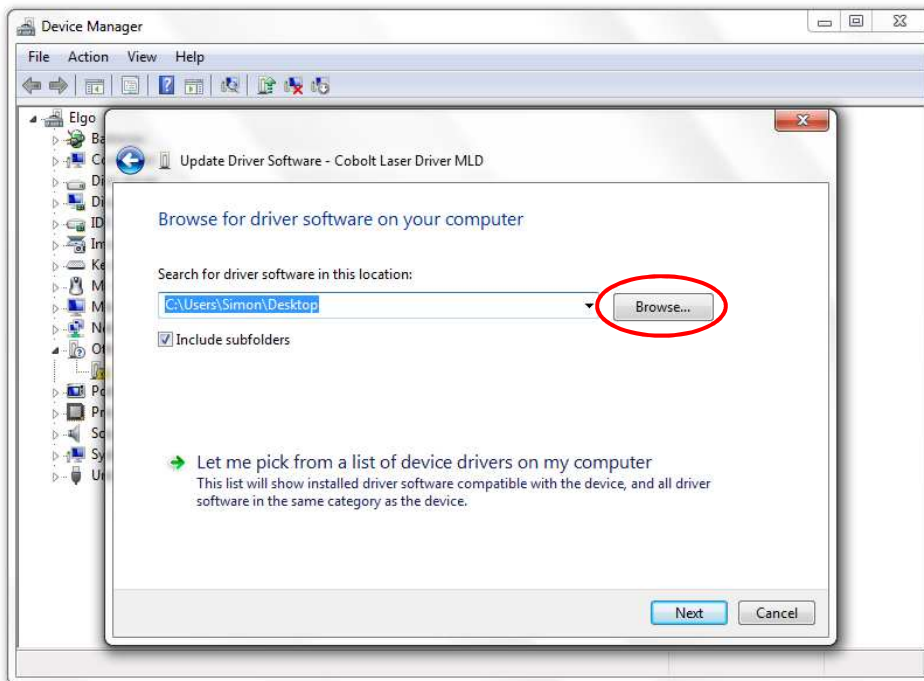

6. Choose the folder and follow the windows instructions to finish the installation.